

## Product Information

### RNase A Solution (10mg/mL)

Catalogue Number	Size
ATR-N707	1 mL (10 mg/mL)

## Product Description

Bovine pancreatic RNase A, a member of the RNase A protein superfamily, is a well-characterized kidney-shaped basic protein consisting of 124 amino acids. In its native state, RNase A forms a homodimer. This endoribonuclease specifically degrades single-stranded RNA at C and U residues by cleaving the phosphodiester bond between the 5'-ribose of a nucleotide and the phosphate group attached to the 3'-ribose of an adjacent pyrimidine nucleotide, resulting in the hydrolysis of the resulting 2', 3'-cyclic phosphate to the corresponding 3'-nucleoside phosphate.

Treatment with RNase A is optional in protocols aiming to remove residual RNA from Genomic DNA Purification Kits to prevent the overestimation of DNA yield due to co-purification of RNA during gDNA extraction. This high-quality enzyme preparation, purified from bovine pancreas, is free of DNase, protease, and DNA contaminants.

## Applications

- **Plasmid and genomic DNA preparation**
- **Removal of RNA from recombinant protein preparations**
- **Ribonuclease protection assays, used in conjunction with RNase T1**
- **Mapping single-base mutations in DNA or RNA**

## Highlights

- **Ready-to-use solution**
- **High-quality enzyme**
- **DNase, protease and DNA-free**

## Inhibition and Inactivation

Inhibitors of ribonucleases include mammalian ribonuclease inhibitor, uridine 2',3'-cyclic vanadate, 5'-diphosphoadenosine 3'-phosphate, 5'-diphosphoadenosine 2'-phosphate, SDS, diethyl pyrocarbonate (DEPC), 4 M guanidinium thiocyanate plus 0.1 M 2-mercaptoethanol, and heavy metal ions.

The most potent inhibitor, mammalian ribonuclease inhibitor, competitively inhibits RNase A. Other inhibitors act through mechanisms such as denaturing DNA, with denatured DNA exhibiting a higher inhibitory effect than native nucleic acids. Additionally, alkylation of His12 or His119, present in the enzyme's active site, can also inhibit RNase A. However, RNase A remains highly active under diverse conditions and is challenging to inactivate. Methods for RNase A inactivation include spin column purification, or phenol/chloroform extraction.

## Note

The recommended concentration range for RNase A is 1 to 100 µg/mL, with the optimal concentration varying based on the specific application. The enzyme exhibits activity across a broad spectrum of reaction conditions. RNase A demonstrates peak activity at a pH range of 7.0 to 7.5. Under low salt concentrations ranging from 0 to 100 mM NaCl, RNase A effectively cleaves both single-stranded and double-stranded RNA, including the RNA strand in RNA-DNA hybrids. However, at NaCl concentrations of 0.3 M or higher, RNase A selectively targets single-stranded RNA for cleavage.

## Storage

Stored at -20°C.

## Shipping

The enzyme is shipped with ice gel.

## Precautions and Disclaimer

This product is intended for research and development (R&D) purposes only and is not suitable for use in drugs, diagnostic procedures, households, or other applications. When handling the product, always wear appropriate laboratory attire, including a lab coat, disposable gloves, and protective

eyewear. When working with radioactive tracers, adhere to standard safety protocols for handling radioactive materials. For detailed safety information, refer to the relevant material safety data sheets (MSDSs), which are available online as PDF files or upon request via email ([info@atrmed.com](mailto:info@atrmed.com)). To the fullest extent permitted by law, ATR-MED Inc. disclaims liability for any special, incidental, indirect, punitive, multiple, or consequential damages arising from or related to the use of this document, including any associated products. By using this product, you acknowledge and agree to all terms and conditions outlined by ATR-MED. All trademarks mentioned herein are the property of ATR-MED unless otherwise specified.

### Limited Product Warranty

ATR-MED® warrants its products in accordance with the terms outlined in ATR-MED's General Terms and Conditions of Sale, available at [www.atrmed.com/terms-and-conditions.html](http://www.atrmed.com/terms-and-conditions.html). For inquiries regarding warranty coverage or any other concerns, please contact ATR-MED via [www.atrmed.com/support](http://www.atrmed.com/support). ATR-MED warrants that, at the time of quality release or subsequent retest date, this product conforms to the information provided in this publication. However, please note that the information contained in this guide is subject to change without prior notice. Purchasers are responsible for determining the suitability of the product(s) for their specific applications. Additional terms and conditions of sale may be found on the invoice or packing slip.

### Trademarks

ATR-MED® is a registered trademark of Acell Teb Rad Med Iranian.